

Operating Instructions Hanging Balance

KERN CH

Version 2.7

07/2009

GB



1 Technical Data

KERN	CH15K20	CH50K50	CH50K100	CH50K500
Readout	20 g	50g	100g	500g
Weighing range	15 kg	50kg	50kg	50kg

2 Fundamental information (general)

2.1 Intended use

The balance you have acquired serves to determine the weighing value of the material to be weighed. It is intended to be used as a "non-automatic" balance, i.e. the material to be weighed is manually and carefully attached at the balance. The weighing value can be read off after a stable weighing value has been obtained.

2.2 Inappropriate use

Do not use the balance for dynamic weighing. In the event that small quantities are removed or added to the material to be weighed, incorrect weighing results can be displayed due to the "stability compensation" in the balance. (Example: Slow draining off of liquid from a container suspended from the balance) Do not attach a continuous load. This can damage the measuring unit as well as the parts, relevant to safety.

Prevent jolts, torsion and oscillation (e.g. by appending slopingly) of all kinds. Be sure to prevent overloading the balance in excess of the stated maximum load (max.), minus any tare weight that may possibly exist. This could damage the balance (risk of breakage).

Important:

- **Always make sure that there are no people or materials below the load that could be injured or damaged!**
- **The balance is not suitable for weighing people. Do not use as baby scales!**
- **The balance does not comply with the medical product law (MPG).**

Never operate the balance in hazardous locations. The series design is not explosion-proof. Structural alterations may not be made to the balance. This can lead to incorrect weighing results, faults concerning safety regulations as well as to destruction of the balance. The balance may only be used in compliance with the described guidelines. Varying areas of application/planned use must be approved by KERN in writing.

2.3 Guarantee

The guarantee is not valid following

- non-observation of our guidelines in the operating instructions
- use outside the described applications
- alteration to or opening of the device
- mechanical damage and damage caused by media, liquids
- natural wear and tear
- inappropriate erection or electric installation
- overloading of the measuring equipment

2.4 Monitoring the test substances

The metrology features of the balance and any possible available adjusting weight must be checked at regular intervals within the scope of quality assurance. For this purpose, the answerable user must define a suitable interval as well as the

nature and scope of this check. Information is available on KERN's home page (www.kern-sohn.com) with regard to the monitoring of balance test substances and the test weights required for this. Test weights and balances can be adjusted quickly and at a reasonable price in KERN's accredited DKD calibration laboratory (return to national normal).

3 Fundamental safety information

- ⚠ Do not use the hanging balance to transport loads.
- ⚠ Prevent jolts, torsion and oscillation (e.g. by appending slopingly) of all kinds.



- ⚠ Never use the hanging balance over the maximum permitted weight (!!Danger of breaking!!).
- ⚠ Always make sure that there are **no living beings or materials** below the load that could be injured or damaged.
- ⚠ The hanging electronic balances from the **KERN CH** series are **only** suitable for hand-held use (see photo). They are **not** suitable for hanging from a mechanical hook, e.g. a crane hook.



3.1 Observe the information in the operating instructions
Please read the operating instructions carefully before erecting and commissioning, even if you already have experience with KERN balances.

3.2 Staff training

The device may only be operated and looked after by trained members of staff.

4 Transport and storage

4.1 Acceptance check

Please check the packaging immediately upon delivery and the device during unpacking for any visible signs of external damage.

4.2 Packaging

Please retain all parts of the original packaging in case it should be necessary to return items at any time. Only the original packaging should be used for return consignments. Before despatch, disconnect all attached cables and loose/movable parts. Apply any intended transport security devices. Secure all parts, e.g. glass windshield, weighing plate, power unit etc., to prevent slipping and damage.

5 Unpacking, installation and commissioning

5.1 Place of installation, place of use

The balance is constructed in such a way that reliable weighing results can be achieved under normal application conditions. By selecting the correct application location for your balance, you will be able to work quickly and precisely.

Therefore please observe the following at the application location:

- Avoid extreme heat as well as temperature fluctuation caused by installing next to a radiator or in the direct sunlight;
- Protect the balance against direct draughts due to open windows and doors;
- Avoid jarring during weighing;
- Protect the balance against high humidity, vapours and dust;
- Do not expose the device to extreme dampness for longer periods of time. Inadmissible bedewing (condensation of air moisture on the device) can occur if a cold device is taken into a significantly warmer environment. In this case, please acclimatise the device for approx. 2 hours at room temperature after it has been disconnected from the batteries.
- Avoid static charging of the material to be weighed, weighing container and windshield.

Major display deviations (incorrect weighing results) are possible if electromagnetic fields occur as well as due to static charging and instable power supply. It is then necessary to change the location.

5.2 Unpacking

Carefully remove the balance from its packaging, remove the plastic wrapping and position the balance in its intended working location.

5.2.1 List of items supplied

Standard accessories:

- Hanging balance
- Battery
- Operating instructions

5.3 Battery operation and change

Push off the battery cover on the back of the balance in the direction of the arrow. Insert 9 V battery. Replace the battery cover. When the batteries are exhausted the message "LO" appears in the balance display. Press the OFF switch and replace the batteries immediately. To preserve battery life the balance switches off automatically after 4 minutes if no weighing has taken place. Deactivate this Auto-OFF function in the menu as follows:#

5.4 The menu

- ⇒ When the balance is switched off, press the **HOLD**-key and keep it pressed
- ⇒ Do not release the **HOLD**-key. Press also the **ON/OFF/TARE**-key and keep it pressed
- ⇒ Keep the **ON/OFF/TARE**-key pressed, however release the **HOLD**-key
- ⇒ Press **HOLD**-key anew
- ⇒ Keep both keys pressed until in the display "tr" appears
- ⇒ Release both keys. The balance is situated in the menu.
- ⇒ Using the ON/OFF/TARE button you can select between the following functions:
 - **tr** (zero tracking) : **ON/OFF**
 - **AF** (Auto off): **ON/OFF**
 - **rST** (Reset to factory setting): **YES/NO**
- ⇒ Confirm with **HOLD**-key
- ⇒ In the display appears the current setting „**ON**“ or „**OFF**“ resp. „**YES**“ or „**NO**“. With help of the **ON/OFF/TARE**-key you can select between „**ON**“ or „**OFF**“ resp. „**YES**“ or „**NO**“. Confirm your selection with the **HOLD**-key. After a short period the balance returns automatically to weighing mode.

6 Adjustment

As the acceleration value due to gravity is not the same at every location on earth, each balance must be coordinated – in compliance with the underlying physical weighing principle – to the existing acceleration due to gravity at its place of location (only if the balance has not already been adjusted to the location in the factory). This adjustment process must be carried out during the initial start-up, after change in location and variation of surrounding temperature. It is also recommendable to adjust the balance periodically during weighing operation in order to obtain exact measured values.

6.1 Adjusting

Using a precision weight, the accuracy of the balance can be checked at any time and adjusted.

Adjustment procedure:

Check that the surrounding conditions are stable. A short warm-up time of about 1 minutes are recommended for stabilisation.

Press "**UNIT**" key approx. 3 seconds, "**CAL**" and then the exact weight of the adjusting weight will appear at the display.

Hang on the adjusting weight on the hook of the hanging scale. "**F**" will appear a short time later, and then turn off.

In case of an adjusting error or a wrong adjusting weight "**E**" appears in the display.

Recommended Adjusting weight: (not included)

CH15K20	CH50K50	CH50K100	CH50K500
10kg (M3)	20kg (M3)	20kg (M3)	20kg (M3)

7 Operation Instruction

7.1 Weighing

- The balance is switched on by pressing the "**ON/OFF/TARE**" key.
- The balance display shows "**0.00**".
- Hang on the weighing object carefully.
- If the weighing object is heavier than the weighing range the display indicates "**E**" (=Overload). Be sure to prevent overloading (chap. 2.2 "Inappropriate use").
- The balance is switched off by pressing the "**ON/OFF/TARE**" key once.

7.2 Weighing units switch-over

Using the "**UNIT**" key you can switch over the display between **N**, **lb** and **kg**. The ► marks the active setting.

7.3 Taring

- Switch the balance on with "**ON/OFF/TARE**" key and wait for "0" to appear in the display. Hang on the tare load carefully and press the "**ON/OFF/TARE**" key again.
- The balance display becomes "**0.00**". The weight of the tare load has however been memorised internally. The item to be weighed can now be placed carefully in the tare container and the weight can be read off.

7.4 HOLD-key / animal weighing / peak value weighing

The HOLD-key may be used to activate various functions. To select and activate these functions, please follow the sequence of operations below:

- Switch on the balance and keep pressing the **HOLD**-key until "H x" appears on the display. The currently activated function is shown above "H1-H6" and by a triangle on the display.
- To select one of the functions, please press the **ON/OFF/TARE** -key. The setting is stored by briefly pressing the **HOLD**-key.

7.4.1 Functions

HOLD function:

The HOLD function **allows** the fixed display of weighing values for longer periods of time.

Depending on the operating range the ON/OFF switch, criteria may be selected from 4 options (H1-H4).

- Selecting the setting "H1" freezes the weighing values for 5 seconds after the HOLD-key was operated.
- Selecting "H2" freezes the weighing values from operating the HOLD-key until another key is pressed.
- Selecting "H3" freezes the weighing values automatically for 5 seconds.
- Selecting "H4" freezes the weighing values automatically until a key is pressed.

HOLD function is set in the following way:

- Switch on the balance by pressing **ON/OFF/TARE** key.
- Press and hold the **HOLD** key until the display shows the required setting (check above „H1–H4“)
- Save the setting by brief pressing the **HOLD** key

Animal weighing function:

This function is suitable for unsteady weighing procedures. A mean value is derived from 16 weighing values.

The animal weighing function is set in the following way:

- Switch on the balance by pressing the **ON/OFF/TARE** key.
- Press and hold the **HOLD** key until „H x“ is shown on the display.
- Press the **ON/OFF/TARE** key repeatedly until „H5“ is shown on the display.
- Save the setting by briefly pressing the **HOLD** key.
- Once again press **HOLD** key.
- On the left side, on top of the display, a triangle will start to blink.
- Hang on the weighing object.
- During the triangle blinking, the weighing values are loaded.
- Then an average weight value is shown consisting of the loaded values.

Peak value function:

This function displays the maximum load value (peak value) of a weighing process.

Measuring frequency: 200ms

Attention:



Under no circumstances may the peak value exceed the stated maximum load of the balance (risk of fracture!!).

How to set the peak value.

- Switch on the balance by pressing the **ON/OFF/TARE** key.
- Press and hold the **HOLD** key until „H x“ is shown on the display.
- Press the **ON/OFF/TARE** repeatedly until „H6“ setting is shown on the display.
- Save the setting by briefly pressing the **HOLD** key.
- Hang on the weighing object.
- The operation of the peak value function is shown by the triangle on top left side of the display.

The negative value (=) is shown right after unloading the balance.

8 Maintenance, upkeep, disposal

8.1 Cleaning

Only use a cloth dampened with mild suds and not aggressive cleaning agents (solvents or similar). Please ensure that fluids are not able to get into the device and rub off using a clean, soft cloth. Loose sample residue/powder can be removed carefully using a brush or hand vacuum cleaner.

8.2 Maintenance, upkeep

The operating company shall ensure that load-bearing components are checked by a competent person at maximum intervals of one year.

These checks can be necessary at intervals shorter than one year depending on the conditions of use. This applies, for example, in the case of particularly frequent use, increased wear and corrosion.

The device may only be opened by trained service engineers authorised by KERN.

Disconnect from the mains supply before opening.

8.3 Disposal

The operating company shall dispose of the packaging and the device in compliance with the valid national or regional law of the operating location.

9 Troubleshooting

The balance should be switched off for a short time following an interruption in the programme sequence. It is then necessary to repeat the weighing process from the beginning.

Interruption	Possible cause
Weight display is not illuminated.	<ul style="list-style-type: none">• The balance is not switched on.• The batteries are wrongly inserted, the batteries are empty• No batteries are attached
The weight display changes continually	<ul style="list-style-type: none">• Draught/air movement• Table/floor vibrations• Vibrations on the balance suspension• Electromagnetic fields / static charging (choose different location/switch off interfering device if possible)
The weighing result is obviously incorrect	<ul style="list-style-type: none">• The balance display is not set to zero• Adjustment is no longer correct.• Great fluctuations in temperature.• Electromagnetic fields / static charging (choose different location/switch off interfering device if possible)

Switch the balance off if other error messages should appear and then switch on again. Contact the manufacturer if the error message does not disappear.